

青霉素-链霉素溶液(100×双抗)

产品简介:

青霉素-链霉素溶液(100×双抗)(Penicillin-Streptomycin Solution, 100×)中含有 10kU/ml 青霉素和 10mg/ml 链霉素, 用 PBS 配制而成, 经过滤除菌, 可直接添加到细胞培养液内用于细胞培养。在细胞培养液中推荐的青霉素的工作浓度为 100U/ml, 链霉素的工作浓度为 0.1mg/ml, 即按照 100 倍稀释本品即可使用; 100ml 本品可配制 10L 细胞培养液。该试剂仅用于科研领域, 不适用于临床诊断或其他用途。

产品组成:

名称 \ 编号	CA0075	Storage
青霉素-链霉素溶液(100×双抗)	100ml	-20℃
使用说明书	1 份	

操作步骤(仅供参考):

参考如下两种方法之一使用:

- 1、在无菌的细胞培养液中直接添加: 按照每 500ml 细胞培养液添加 5ml 的比例加入青霉素-链霉素溶液(100×双抗), 混匀即可使用。
- 2、配制细胞培养液时加入, 然后再过滤除菌: 按照每配制 1L 细胞培养液加入 10ml 的比例加入本品, 配制完成后过滤除菌即可使用。

注意事项:

- 1、尽量减少反复冻融的次数, 以免效率下降。
- 2、试剂开封后请尽快使用, 以防影响后续实验效果。
- 3、注意无菌操作, 尽量避免污染。
- 4、为了您的安全和健康, 请穿实验服并戴一次性手套操作。

有效期: 12 个月有效。低温运输, -20℃保存。

相关产品:

产品编号	产品名称
CA0005	氨苄青霉素溶液(Ampicillin,50mg/ml)
CC0005	1×PBS 缓冲液(无钙镁)
CC0122	无酶细胞消化液
CC0128	胰蛋白酶-EDTA 溶液(0.25%:0.02%,含酚红)
CT0022	Cell Counting Kit-8 (CCK-8 试剂盒)
CT0027	乳酸脱氢酶(LDH)细胞毒性检测试剂盒
CZ0016	LipoLeagene2000 转染试剂
DA0010	Hoechst33258 染色液
DA0039	AO/EB 双荧光染色试剂盒
DF0135	组织细胞固定液(4% PFA)

文献引用:

- 1、Yudi Deng,Zhijie Jiang,Yincheng Jin,et al.Reinforcing vascular normalization therapy with a bi-directional nano-system to achieve therapeutic-friendly tumor microenvironment.JOURNAL OF CONTROLLED RELEASE E.October 2021.10.1016/j.jconrel.2021.10.016.(IF 9.776)
 - 2、Wu Peipei,Tang Yuting,Jin Can,et al.Neutrophil membrane engineered HucMSC sEVs alleviate cisplatin-induced AKI by enhancing cellular uptake and targeting.JOURNAL OF NANOBIO TECHNOLOGY.August 2022. 10.1186/s12951-022January574-8.(IF 9.429)
 - 3、Xuechen Zhu,Pan Wang,Jiale Wei,et al.Lysosomal degradation of the maternal dorsal determinant Hwa safeguards dorsal body axis formation.EMBO REPORTS.October 2021.10.15252/embr.202153185.(IF 8.807)
 - 4、Huang Yusheng,Xia Lei,Tan Xiangwu,et al.Molecular mechanism of lncRNA SNHG12 in immune escape of non-small cell lung cancer through the HuR/PD-L1/USP8 axis.CELLULAR & MOLECULAR BIOLOGY LETTERS.June 2022.10.1186/s11658-022-00343-7.(IF 8.702)
 - 5、Feng Ouyang,Yu Liu,Li Zhang,et al.Hyaluronic acid-targeted dual-bubble/photothermal-driven nanomissile for enhanced “four-in-one” anti-tumor strategy.INTERNATIONAL JOURNAL OF BIOLOGICAL MACROMOLECULES.July 2024.10.1016/j.ijbiomac.2024.133997.(IF 7.7)
 - 6、Qiuhong Xiong,Huimin Sun,Wenxiu Xing,et al.WDR45 mutation dysregulates iron homeostasis by promoting the chaperone-mediated autophagic degradation of ferritin heavy chain in an ER stress/p38 dependent mechanism.FREE RADICAL BIOLOGY AND MEDICINE.March 2023.10.1016/j.freeradbiomed.2023.03.012.(IF 7.4)
 - 7、F. Ouyang,L. Zhang,Y. Liu,et al.Environmentally responsive multistage gas/enzyme/NIR-driven chemotactic transport nanorockets for single wavelength procedural combination therapy.Materials Today Chemistry.April 2023.10.1016/j.mtchem.2023.101533.(IF 7.3)
 - 8、Yang Yakun,Qi Jiaying,Zhang Muqing,et al.The cardioprotective effects and mechanisms of naringenin in myocardial ischemia based on network pharmacology and experiment verification.Frontiers in Pharmacology.September 2022.10.3389/fphar.2022.954555.(IF 5.988)
 - 9、Guo Haohao,Ma Runsheng,Zhang Yifei,et al.Ibuprofen inhibits anaplastic thyroid cells in vivo and in vitro by triggering NLRP3-ASC-GSDMD-dependent pyroptosis.INFLAMMOPHARMACOLOGY.November 2023.10.1007/s10787-023January379-7.(IF 5.8)
 - 10、Zixin Deng,Defu Liu,Yi Xiong,et al.Preparation of a hydroxyapatite-silver gradient bioactive ceramic coating with porous structure by laser cladding: A study of in vitro bioactivity.CERAMICS INTERNATIONAL.July 2022.10.1016/j.ceramint.2022.06.327.(IF 5.532)
 - 11、Tianshu Xu,Ping He,So namWangdu,et al.Revealing the improvement of diabetes by Si Wei Jiang Huang Tang San through ERK/HIF1 α signaling pathway via network pharmacology.JOURNAL OF ETHNOPHARMACOLOGY.September 2023.10.1016/j.jep.2023.117254.(IF 5.4)
 - 12、Fen Wei,Xiaxi Zhang,Ping Cui,et al.Cell-based 3D bionic screening by mimicking the drug-receptor interaction environment in vivo.Journal of Materials Chemistry B.December 2020.10.1039/D0TB02661A.(IF 5.344)
 - 13、Lijun Zhang,Zhenning Dai,Shanshan Shi,et al.SIRT3 and SIRT4 double-genes remodeled the mitochondrial network to induce hepatocellular carcinoma cell line differentiation and suppress malignant phenotypes.BIOCHEMICAL PHARMACOLOGY.March 2024.10.1016/j.bcp.2024.116168.(IF 5.3)
 - 14、Yakun Yang,jiaying Qi,Hua Wang,et al.Ameliorative effects of Danshensu from the functional food Salvia miltiorrhiza against arsenic trioxide-induced cardiac toxicity in vivo and in vitro: Involvement of inhibiting the AKT/IKK/NF- κ B signaling pathway.Journal of Functional Foods.October 2022.10.1016/j.jff.2022.105304.(IF 5.223)
 - 15、Jinjing Xia,Hao Bai,Bo Yan,et al.Mimicking the BIM BH3 domain overcomes resistance to EGFR tyrosine kinase inhibitors in EGFR-mutant non-small cell lung cancer.Oncotarget.July 2017.10.18632/oncotarget.19411.(IF 5.168)
 - 16、Yan Zhou,Lu-Lu Wen,Yan-Fei Li,et al.Exosomes derived from bone marrow mesenchymal stem cells protect the injured spinal cord by inhibiting pericyte pyroptosis.Neural Regeneration Research.June 2021.10.4103/1673-5374.314323.(IF 5.135)
 - 17、Xue Fumin,An Chao,Chen Lixiang,et al.4.1B suppresses cancer cell proliferation by binding to EGFR P13 region of intracellular juxtamembrane segment.Cell Communication and Signaling.September 2019.10.1186/s12964January9-0431-6.(IF 5.111)
 - 18、Deqian Kong,Yue Liu,Li Li,et al.Astaxanthin ameliorates oxidative stress in lens epithelial cells by regulating GPX4 and ferroptosis.CHEMICO-BIOLOGICAL INTERACTIONS.August 2023.10.1016/j.cbi.2023.110684.(IF 5.1)
-

- 19、Ying Chen,Bolin Wu,Yameng Hao,et al.Structure-activity relationship studies of (E)-3,4-dihydroxystyryl alkyl sulfones as novel neuroprotective agents based on improved antioxidant, anti-inflammatory activities and BBB permeability.EUROPEAN JOURNAL OF MEDICINAL CHEMISTRY.March 2019.10.1016/j.ejmech.2019.03.044.(IF 4.833)
- 20、Jiaying Qi,Haoying Li,Yakun Yang,et al.Mechanistic insights into the ameliorative effects of hypoxia-induced myocardial injury by Corydalis yanhusuo total alkaloids: based on network pharmacology and experiment verification.Frontiers in Pharmacology.January 2024.10.3389/fphar.2023.1275558.(IF 4.4)
- 21、Xiong Qihong,Sun Huimin,Wang Yanlin,et al.Lipid droplet accumulation in Wdr45-deficient cells caused by impairment of chaperone-mediated autophagic degradation of Fasn.Lipids in Health and Disease.March 2024.10.1186/s12944-024-02088-y.(IF 3.9)
- 22、Shumiao Zhang,Fan Yang,Rui Shi,et al.Activation of κ -opioid receptor inhibits inflammatory response induced by sodium palmitate in human umbilical vein endothelial cells.CYTOKINE.July 2021.10.1016/j.cyto.2021.155659.(IF 3.861)
- 23、Feng Ouyang,Li Zhao,et al.Photothermally driven degradable nanomissile for single-wavelength phototherapy and hypoxia-activated chemotherapy.INORGANIC CHEMISTRY COMMUNICATIONS.August 2023.10.1016/j.inoche.2023.111310.(IF 3.8)
- 24、Yun-xia An,Yi-jun Shang,Zhi-wei Xu,et al.STAT3-induced long noncoding RNA LINC00668 promotes migration and invasion of non-small cell lung cancer via the miR-193a/KLF7 axis.BIOMEDICINE & PHARMACOTHERAPY.May 2019.10.1016/j.biopha.2019.109023.(IF 3.743)
- 25、Han Biqu,Lv Zhanjun,Han Xuemin,et al.Harmful Effects of Inorganic Mercury Exposure on Kidney Cells: Mitochondrial Dynamics Disorder and Excessive Oxidative Stress.BIOLOGICAL TRACE ELEMENT RESEARCH.May 2021.10.1007/s12011-021-02766-3.(IF 3.738)
- 26、Dan-Dan Sun,Xue Wu,Shi-Chen Lin,et al.Anti-apoptosis and anti-inflammation activity of circ_0097010 downregulation in lipopolysaccharide-stimulated periodontal ligament cells by miR-769-5p/Krüppel like factor 6 axis.Journal of Dental Sciences.June 2022.10.1016/j.jds.2022.04.024.(IF 3.719)
- 27、Hairui Zhou,Jingru Wu,Yu Gong,et al.Isoquercetin alleviates sleep deprivation dependent hippocampal neurons damage by suppressing NLRP3-induced pyroptosis.IMMUNOPHARMACOLOGY AND IMMUNOTOXICOLOGY.June 2022.10.1080/08923973.2022.2082976.(IF 3.712)
- 28、Song Gao,Gao, Jianchao Wang, Wang, Shujian Tian,et al.miR-9 depletion suppresses the proliferation of osteosarcoma cells by targeting p16.INTERNATIONAL JOURNAL OF ONCOLOGY.April 2019.10.3892/ijo.2019.4783.(IF 3.571)
- 29、Yun Huang,Qing Peng,Hai-Yan Li,et al.Novel sericin-based hepatocyte serum-free medium and sericin's effect on hepatocyte transcriptome.WORLD JOURNAL OF GASTROENTEROLOGY.August 2018.10.3748/wjg.v24.i30.3398.(IF 3.3)
- 30、Ming Fang,Yanfei Li,Yingbiao Wu,et al.miR-185 silencing promotes the progression of atherosclerosis via targeting stromal interaction molecule 1.CELL CYCLE.March 2019.10.1080/15384101.2019.1580493.(IF 3.259)
- 31、Haohao Guo,Ziyang Wang,Keyu Yin,et al.Sciellin promotes the development and progression of thyroid cancer through the JAK2/STAT3 signaling pathway.MOLECULAR CARCINOGENESIS.February 2024.10.1002/mc.23682.(IF 3)
- 32、Feng Qu,Xuezheng Shen,Ketao Wang,et al.Tenogenic differentiation of human tendon-derived stem cells induced by long non-coding RNA LINCMD1 via miR-342-3p/EGR1 axis.CONNECTIVE TISSUE RESEARCH.June 2023.10.1080/03008207.2023.2217258.(IF 2.9)
- 33、Guangyu Sun,Yanmin Song,Congxian Li,et al.MTCH2 promotes the malignant progression of ovarian cancer through the upregulation of AIMP2 expression levels, mitochondrial dysfunction and by mediating energy metabolism.Oncology Letters.August 2024.10.3892/ol.2024.14625.(IF 2.5)
- 34、Yan-Ci Li,Ke Tian,Li-Juan Sun,et al.A new hexacyclic triterpene acid from the roots of Euscaphis japonica and its inhibitory activity on triglyceride accumulation.FITOTERAPIA.January 2016.10.1016/j.fitote.2016.01.016.(IF 2.408)
- 35、Yujing Fan,et al.miR-205 suppresses cell migration, invasion and EMT of colon cancer by targeting mouse double minute 4.Molecular Medicine Reports.May 2020.10.3892/mmr.2020.11150.(IF 2.1)
- 36、Peihong Shen,Shen, Junfeng Chen, Chen,et al.The protective effects of total paeony glycoside on ischemia/reperfusion injury in H9C2 cells via inhibition of the PI3K/Akt signaling pathway.Molecular Medicine Reports.July 2018.10.3892/mmr.2018.9335.(IF 1.922)
- 37、Chunxia Yu,Weimin Wang,et al.Hirudin Protects Ang II-Induced Myocardial Fibroblasts Fibrosis by Inhibiting the Extracellular Signal-Regulated Kinase1/2 (ERK1/2) Pathway.MEDICAL SCIENCE MONITOR.Septemb

-
- er 2018.10.12659/MSM.909044.(IF 1.894)
- 38、 Shushu Wang,Yueyang Liang,Wenxiao Chang,et al.Triple Negative Breast Cancer Depends on Sphingosine Kinase 1 (SphK1)/Sphingosine-1-Phosphate (S1P)/Sphingosine 1-Phosphate Receptor 3 (S1PR3)/Notch Signaling for Metastasis.MEDICAL SCIENCE MONITOR.April 2018.10.12659/MSM.905833.(IF 1.894)
- 39、 Xuebing Wang,Xiao Li,Xuefang Wang,et al.Experimental study of Forsythoside A on prevention and treatment of avian infectious bronchitis.RESEARCH IN VETERINARY SCIENCE.November 2020.10.1016/j.rvsc.2020.11.009.(IF 1.892)
- 40、 Jun Zhang,Zhang, Rongming Xu, Xu, Lei Wu,et al.Expression and function of Toll- like receptors in peripheral blood mononuclear cells in patients with ankylosing spondylitis.Molecular Medicine Reports.September 2019.10.3892/mmr.2019.10631.(IF 1.851)
- 41、 Liu Yayun,Xi Yizhe,Chen Gang,et al.URG4 mediates cell proliferation and cell cycle in osteosarcoma via GSK3 β / β -catenin/cyclin D1 signaling pathway.Journal of Orthopaedic Surgery and Research.June 2020.10.1186/s13018-020January681-y.(IF 1.777)
- 42、 Xuan Kan,Yanan Sun,Jianguang Lu,et al.Co- inhibition of miRNA- 21 and miRNA- 221 induces apoptosis by enhancing the p53- mediated expression of pro- apoptotic miRNAs in laryngeal squamous cell carcinoma.Molecular Medicine Reports.March 2016.10.3892/mmr.2016.5048.(IF 1.559)
- 43、 Yanyan Wang,Qiang Guo,Zhongbin Cheng,et al.New saikosaponins from the roots of Bupleurum chinense. Phytochemistry Letters.July 2017.10.1016/j.phytol.2017.06.005.(IF 1.418)
- 44、 Huaiyu Tong,Yan Wang,Xuechun Lu,et al.On the preparation of transferrin modified artesunate nanoliposomes and their glioma-targeting treatment in-vitro and in-vivo.International Journal of Clinical and Experimental Medicine.December 2015.PMID:26885179.(IF 1.277)

注：更多使用本产品的文献请参考产品网页